

for  
**10mL**  
5mL  
pipette


**No electric power  
required at all!**

**Repetitive Pipette**


# Peggy



# Aspirate the sample all at once, dispensing its necessary volume repetitively.

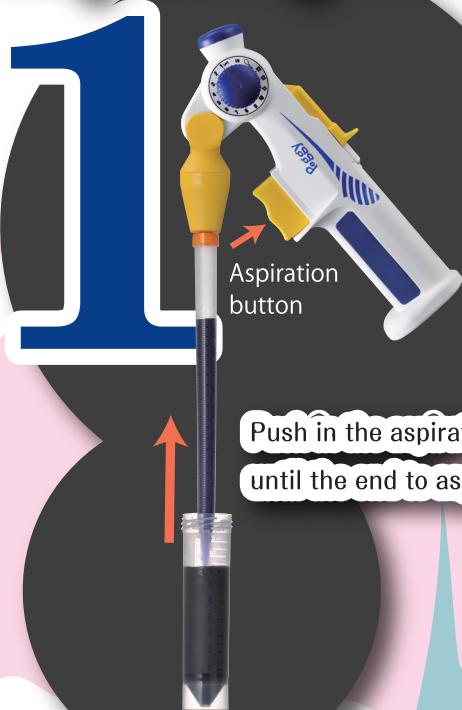


Push in the aspiration button once slowly until the end to aspirate the set volume.

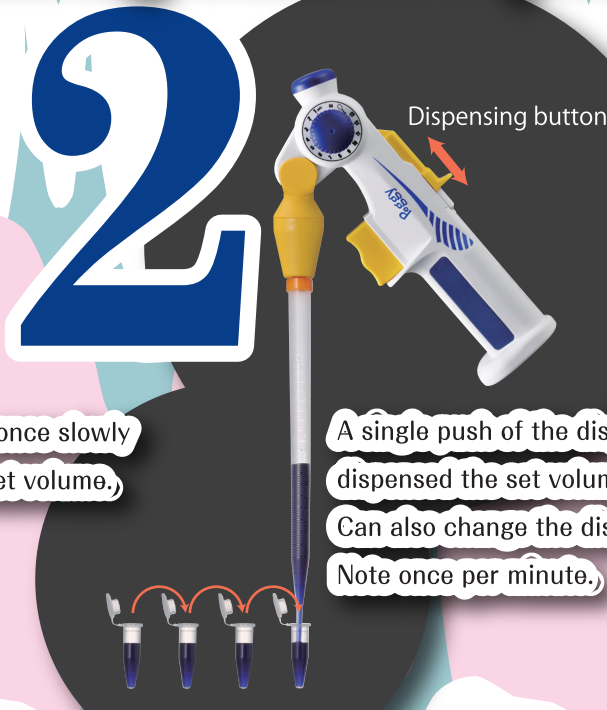

A single push of the dispensing button dispensed the set volume.  
Can also change the dispensing capacity Note once per minute.


## Setting aspiration volume

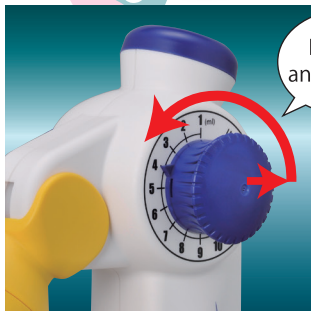

Aspiration can be set in every 1mL for setting range 1-13mL.


Note: 13mL can be dispensed in case that Watson 10mL Serological Pipette is used.

## Setting dispensing volume

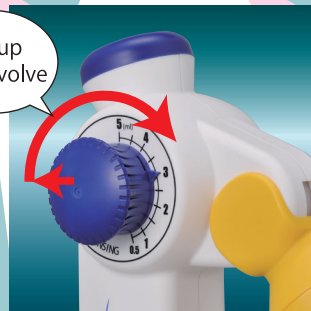


Dispensing can be set in every 0.5mL for setting range 0.5-5mL.

## Flexible nozzle-angle

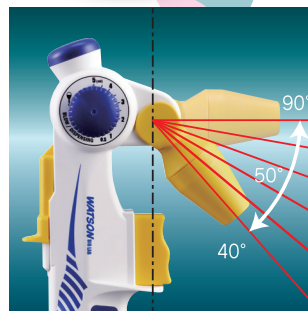

Adjustable nozzle angle in every 10 degrees up 50 degrees.

## Simple maintenance

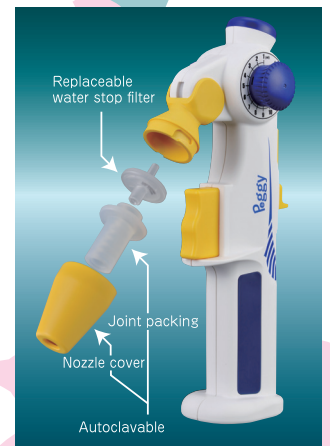

Easy to remove & set spare parts  
- Nozzle cover: Autoclavable  
- Joint packing: Autoclavable  
- water stop filter: Replaceable

### Precaution in case of using 5mL serological pipette:

Please DO NOT set aspiration volume exceeding 5mL in case that 5mL serological pipette is used. It may cause an accidental liquid suction leading to malfunction on Peggy.

Cat.No.	Item	Aspiration Volume		Discharge Volume		Packaging
		max.	min.	max.	min.	
PG-01	PEGGY-01 Repetitive Pipette "Peggy"	13 mL	1 mL	5 mL	0.5 mL	1 unit / box

Something Different.

**WATSON** BIO LAB  
MADE IN JAPAN SINCE 1988

Manufactured and sold by:

**FUKAEKASEI** CO., LTD.

Head Office: 2-2-7 Murotani, Nishi-ku, Kobe  
651-2241, JAPAN  
TEL +81-78-991-4477  
FAX +81-78-991-4491

E-mail: [info@watson.co.jp](mailto:info@watson.co.jp)

<https://watsonbiolab.com>

※The specification and the price etc. of the product might change without a previous notice.